How to access data generated on the Accela LC - LTQ system and to obtain the required software

Connect to UCL Virtual Private Network (VPN)

https://www.ucl.ac.uk/isd/services/get-connected/ucl-virtual-private-network-vpn

Go to windows

Tools

Select map network drive

💮 🤏 Map N	Jetwork Drive
What n Specify th	etwork folder would you like to map? e drive letter for the connection and the folder that you want to connect to:
Drive: Folder:	Y: ■ Browse Example: \\server\share Image: Reconnect at logon Image: Connect using different credentials Connect to a Web site that you can use to store your documents and pictures.
	Finish Cancel

Drive does not matter

Type in the Folder <u>\\storage.chem.ucl.ac.uk\ms</u> or <u>\\128.40.177.194\ms</u>

Click tick on Reconnect at logon

Click tick on Connect using different credentials

It will ask user name and password

Type under user name: ms-ro

Type under password: frollyug

You will see the following window

Eile Home Share	View				- a x
Pin to Quick Copy Paste	6 Cut Copy path Paste shortcut Paste shortcut	New item •	Properties	Select all Select none Invert selection	
Clipboard	Organise	New	Open	Select	
🗹 📙 🔻					
← → < ↑ 💄 > Ne	etwork > storage.chem.ucl.ac.uk > ms >				✓ Ů Search ms P
Windows (C:)	Name ^	Date modified	Туре	Size	Thermo GC-MS
TOCHIDA EVE (C.	backup	17/12/2017 08:18	File folder		File folder
COMBA EXT (E	LCT Premier XE	16/03/2020 15:11	File folder		
GC-MS Archive E	MassHunterData	07/01/2020 13:19	File folder		
C-MS Archive Dr	AT900	08/08/2018 12:14	File folder		
📕 data	Orbitrap	20/05/2020 11:24	File folder		
Seagate	pub	10/03/2020 10:39	File folder		Data and Jillington DE (01/2010 45-40
UserGuides	Thormo Accolate ITO	20/01/2018 09:20	File folder		Date modified: 23/01/2018 13:49
	Thermo Accelato-LTQ	29/01/2020 14:08	File folder		
IOSHIBA EXT (E:)	TOEChecktune Positive 20191120 104	20/11/2019 11:00	File folder		
2018		19/12/2017 14:21	File folder		
Agilent MassHur	Waters Acquity UPLC-SOD	17/02/2020 14:57	File folder		
FORMS	Waters MALDI-TOF	17/01/2020 14:06	File folder		
how to book inst	D.pdf	12/07/2019 16:20	PDF File	627 KB	
March 2020	grml64-full_2018.12.iso	22/05/2020 23:09	Disc Image File	688,384 KB	
Mass Spectrome	TOFAutotune_Positive_20191021_1627	20/11/2019 11:58	Compressed (zipp	22 KB	
MassHunter	TOFChecktune_Positive_20191120_104	20/11/2019 11:07	Compressed (zipp	23 KB	
Master project 2					
Merel_2019					
Orbitrap results					
Softwares					
Techniques					
Xcalibur					
I Network					
17 items 1 item selected					
		—	<u> </u>		
Type here	to search O		C 🔹 🗖	<u>a</u> 💿	C V 🗤 💴 V V V V V V V V V V V V V V V V V V

Your data will be in the folder Thermo Accela LC-LTQ and stored by year and month when the samples were analysed

Eile Home Share	View						- ¤ ×
Pin to Quick Copy Paste	Cut Copy path Paste shortcut Delete Rename	New item •	Properties	Select all Select none Invert selection			
Clipboard	Organise	New	Open	Select			
🖌 📔 🔺							
← → ヾ ↑ 🚑 > N	etwork > storage.chem.ucl.ac.uk > ms >						✓ ひ Search ms ル
🐛 Windows (C:) 🔷	Name ^	Date modified	Туре	Size			Thermo AccelaLC-LTQ
RECOVERY (D:)	backup	17/12/2017 08:18	File folder				File folder
🥪 TOSHIBA EXT (E:	LCT Premier XE	16/03/2020 15:11	File folder				
GC-MS Archive [MassHunterData	07/01/2020 13:19	File folder				The Rest of States of States
CC MS Archive Dr	AT900	08/08/2018 12:14	File folder				T 1
data	Orbitrap	20/05/2020 11:24	File folder				
Gata	📜 pub	10/03/2020 10:39	File folder				and the state of t
seagate	software	18/01/2018 09:26	File folder				Date modified: 29/01/2020 14:08
UserGuides	🗹 📜 Thermo AccelaLC-LTQ	29/01/2020 14:08	File folder				
🧼 TOSHIBA EXT (E:)	Thermo GC-MS	25/01/2018 15:49	File folder				
2018	TOFChecktune_Positive_20191120_104	20/11/2019 11:00	File folder				
Agilent MassHur	tools	19/12/2017 14:21	File folder				
FORMS	Waters Acquity UPLC-SQD	17/02/2020 14:57	File folder				
how to book inst	Waters MALDI-TOF	17/01/2020 14:06	File folder				
March 2020	D.pdf	12/07/2019 16:20	PDF File	627 KB			
Mass Spectrome	Grmi64-101_2018.12.IS0 TOFAutoture Peritius 20101021 1627	22/05/2020 23:09	Compressed (zing	000,304 ND			
MassHunter	TOPAdotale_rositive_20191021_1021	20/11/2019 11:08	Compressed (zipp	22 KB			
Master project 2		20/11/2015 11:07	compressed (cip)	- 23 KU			
Merel 2019							
Orbitran results							
Contrap results							
Softwares							
Techniques							
Xcalibur							
🕩 Network							
×							
17 items 1 item selected	1						822
1 O Turcher	to count	5	o -	2 📼	6 6 *	एम द्यम	
> Type here	to search O	<u> </u>	<u> </u>		<u> </u>		😗 ^ 😑 🚰 🖼 03/06/2020 📑

Pie to Guide. Copy Paste shortcat More Copy Paste shortcat More Copy Paste shortcat Pie to Guide. Copy paste Pie to Guide. Copy paste Pie to Guide. Copy Pie to Guide. Copy		
Clipboard Organise New Open Select		
🗹 📜 🔻		
← → v ↑ 📕 > Network > storage.chem.ucl.ac.uk > ms > Thermo AccelaLC-LTQ. v U	Search T	h_ P
L Windows (C) ^ _ Name _ Date modified Type Size		
RECOVERY (D:) 2017 from Aug 01/11/2017 11:22 File folder		
TOSHIBA EXT (E: 2019 15/11/2019 13:48 File folder		
C-MS Archive E 2020 05/03/2020 10:14 File folder		
C GC-MS Archive Dr 😥 MSMS library Eloippt 09/03/2018 10.52 Microsoft PowerPo 278 KB		
data		
Seagate		
UserGuides		
TOSHIBA EXT (E)		
2018		
Aglient MassHur		
FORMS		
how to book inst		
March 2020		
Mass Spectrome		
MassHunter		
Kaster project 2		
. Merel 2019		
Orbitrap results		
Softwares		
Cechniques		
Zcalibur		
😻 Network		
4 items		800 K
🖽 🔎 Type here to search O 😰 🚖 🤮 🏦 🔝 🕘 🚳 🖉 🤌 🔹 🕅 💷 🎯 🖓	12:41 03/06/2020	5

Download Xcalibur software for data analysis

Click software (this software runs on Windows XP but should work also on Windows 7, 8, 9, 10, if you have any problems pelase let me know)

ms					- ¤ ×
Pin to Quick Copy Paste	View 6 Cut Copy path Paste shortcut Move Copy to to t	New item •	Properties	Select all	A
Clipboard	Organise	New	Open	Select	
🖌 📜 🔺					
← → ~ ↑ 🕹 > Ne	etwork > storage.chem.ucl.ac.uk > ms >				✓ U Search ms P
😓 Windows (C:) 🔷	Name ^	Date modified	Туре	Size	software
RECOVERY (D:)	backup	17/12/2017 09-19	File folder		File folder
TOSHIBA EXT (E:	ICT Premier YE	16/02/2020 15:11	File folder		
GC-MS Archive E	MassHunterData	07/01/2020 13:19	File folder		
	MAT900	08/08/2018 12:14	File folder		
GC-MS Archive Dr	Orbitrap	20/05/2020 11:24	File folder		
📕 data	pub	10/03/2020 10:39	File folder		
Seagate	Software	18/01/2018 09:26	File folder		Date modified: 18/01/2018 09:26
UserGuides	🗌 📕 Thermo AccelaLC-LTQ	29/01/2020 14:08	File folder		
TOSHIBA EXT (E:)	Thermo GC-MS	25/01/2018 15:49	File folder		
2018	TOFChecktune_Positive_20191120_104	20/11/2019 11:00	File folder		
Agilent MassHur	tools	19/12/2017 14:21	File folder		
	Waters Acquity UPLC-SQD	17/02/2020 14:57	File folder		
have been ind	Waters MALDI-TOF	17/01/2020 14:06	File folder		
Now to book insi	D.pdf	12/07/2019 16:20	PDF File	627 KB	
March 2020	grml64-full_2018.12.iso	22/05/2020 23:09	Disc Image File	688,384 KB	
Mass Spectrome	TOFAutotune_Positive_20191021_1627	20/11/2019 11:58	Compressed (zip)	p 22 KB	
MassHunter	TOFChecktune_Positive_20191120_104	20/11/2019 11:07	Compressed (zip)	р 23 КВ	
Master project 2					
Merel_2019					
Orbitrap results					
Softwares					
Techniques					
Xcalibur					
Network					Ahern, mentioned General in General E Student Health and Wellbeing Community
17 items 1 item selected	1				
		-	· · · ·		- 1529 -

Click software and the Xcalibur

Then click Foundation and run ThermoPlatform

Foundation File Home Share	View								- 🗆 X ^ 🕑
Pin to Quick Copy Paste	Cut Copy path Paste shortcut	Move Copy to * to *	Mew item *	Properties	Select all Select none Invert selection				
Clipboard		Organise	New	Open	Select				
🗹 📜 📼									
← → × ↑ 📜 > Ne	twork ≥ stora	ge.chem.ucl.ac.uk > ms > s	oftware > Xcalibur > Found	ation >					V 🖸 Search Fo 🔎
Windows (C:)	Name	^	Date modified	Туре	Size			2 items	
 RECOVERY (D:) TOSHIBA EXT (E: GC-MS Archive E 	📕 Man 🛃 Ther	uals moPlatform.exe	11/01/2018 09:41 16/08/2011 22:45	File folder Application	15,118 KB				
GC-MS Archive Dr data Seagate									L
TOSHIBA EXT (E:) 2018 Agilent MassHur FORMS how to book inst Massh 2020									
Mass Spectrome MassFlunter Master project 2 Merel_2019									
Softwares Techniques Xcalibur									
Network 2 items									(iii) 🛤
E $\label{eq:powerserver}$ Type here	to search		0 😰 💼	e 💼 토	i a 💿	<i>e</i> 💿	🐞 🚺 🚺	(2) ^ =	a 🖘 🎞 15:31

After click Xcalibur and run XcaliburSetup

Xcalibur File Home Share Pin to Quick Copy Paste Clipboard	View Cut Copy path Paste shortcut	More Copy to to Organise	ne New item * Dessy access * Folder New	Properties Open	Select all Select none Invert selection Select					- 0 X
$\leftarrow \rightarrow \lor \uparrow \blacksquare > Ne$	etwork > storag	ge.chem.ucl.ac.uk > ms > so	ftware > Xcalibur > Xcalib	ur						Search Xc P
Uindows (C:)	Name	^	Date modified	Туре	Size				2 items	
RECOVERY (D:)	📜 Manu	als	11/01/2018 09:40	File folder						
TOSHIBA EXT (E: GC-MS Archive E GC-MS Archive Dr	🛃 Xcalit	burSetup.exe	12/08/2011 19:22	Application	57,640 KB					
data										
Seagate										
UserGuides										
TOSHIBA EXT (E:)										
2018										
Agilent MassHur										
FORMS										
how to book inst										
March 2020										
Mass Spectrome										
MassHunter										
Master project 2										
Merel_2019										
Orbitrap results										
Softwares										
Techniques										
Xcalibur										
Intwork										
~										
2 items										
H \show Type here	to search		0 😰 💼	e 💼 🔚	I a 🗠	🩋 📀	16	×1	🕜 ^ 👄 🍝 😁	15:32 01/06/2020

If you have any questions please email to the mass spec staff