# How to access your data generated on the open access MALDI-ToF mass spectrometer

- Connect to UCL Virtual Private Network (VPN)
- https://www.ucl.ac.uk/isd/services/get-connected/ucl-virtual-private-network-vpn
- Go to windows
- Tools
- Select map network drive

🕞 🤏 Map N	etwork Drive
What ne Specify the	etwork folder would you like to map?
Drive: Folder:	Y: Image: Browse   Example: \\server\share Browse   Reconnect at logon Connect using different credentials   Connect to a Web site that you can use to store your documents and pictures.
	Finish Cancel

- Drive does not matter
- Type in the Folder <u>\\storage.chem.ucl.ac.uk\ms</u> or <u>\\128.40.177.194\ms</u>
- Click tick on Reconnect at logon
- Click tick on Connect using different credentials
- It will ask user name and password
- Type under user name: ms-ro

- Type under password: frollyug
- You will see the following window



- Click on Waters MALDI-TOF
- DATA saved by date/year when MALDI analyses were done

#### For example:-

Data 2016	View						- a x
Pin to Quick Copy Paste	Cut Copy path Paste shortcut Delete Renz	me New folder	Properties	Select all Select none			
Clipboard	Organise	New	Open	Select			
🗹 📜 🔻							
← → × ↑ 📕 > N	etwork > storage.chem.ucl.ac.uk > ms > V	Vaters MALDI-TOF > Data 20	16				V Search Da A
🐛 Windows (C:) 🔷	Name ^	Date modified	Туре	Size			January 2016
RECOVERY (D:)	April 2016	28/04/2016 11:18	File folder				File folder
🧼 TOSHIBA EXT (E:	August 2016	26/08/2016 16:24	File folder				
GC-MS Archive E	December 2016	20/12/2016 16:50	File folder				
In columnities De	February 2016	26/02/2016 15:39	File folder				
GC-MS Archive Dr	January 2016	28/01/2016 15:45	File folder				
data	July 2016	27/07/2016 16:33	File folder				
Seagate	🧯 June 2016	07/06/2016 15:57	File folder				Date modified: 28/01/2016 15:45
UserGuides	March 2016	24/03/2016 11:02	File folder				
TOSHIBA EXT (E:)	May 2016	27/05/2016 10:57	File folder				
2018	November 2016	29/11/2016 14:27	File folder				
Agilent MassHur	October 2016	28/10/2016 11:25	File folder				
FORMS	September 2016	05/10/2016 14:27	File folder				
how to book inst							
March 2020							
Mass Spectrome							
Marshunter							
Master project 2							
Marel 2010							
Orbitrap results							
Cathurana							
Sottwares							
Techniques							
Xcalibur							
🥩 Network							
~							
12 items 1 item selected	d						(800 F
+ P Type here	to search	0 😰 🚖	2 🔒 📮	a 👼	A 💿 💏	x x	🙆 ^ 🛥 🖕 🖓 🛄 🛃

#### Download MassLynx software for data analysis

Click software (this software runs on Windows XP but should work also on Windows 7, 8, 9, 10, if you have any problems pelase let me know)

									X
Compute	er ▶ ms (\\128.40.177.194) (Z:) ▶ software ▶					Search software	_	_	Q
Grganize - Burn	s Help						8== 🗸	FI	
A	Name	Date modified	Type	Size					•
Pavorites	A silest Massiliustes Bis Casting B 07.00	04/01/2019 16-54	Els falsia						
Develop	Agilent MassHunter BioConfirm B.07.00	04/01/2018 10:54	File folder						
Recent Places	Agilent MassHunter Qualitative Analysis	20/12/2017 10:55	File folder						
A Recent Fraces	Massiver 41	02/01/2018 14:02	File folder						
🕞 Libraries	MassLynx 4.1	02/01/2010 14:02	The folder						
Documents									
A Music									
Pictures									
Videos									
Computer									
A 05 (C)									
TOSHIBA EXT (E:)									
TOSHIBA (F:)									
🖵 ms (\\128.40.177.19									
🙀 Network									

Click MassLynx folder and open CD1

Select setup.exe and allow to run this set-up

Click NEXT on each window appeared and accept the virus check and accept the licence question

	Description of the description of the local		And in case of the local division of the loc					
🕞 🗢 📙 🕨 Compu	uter • ms (\\128.40.177.194) (Z:) • software	MassLynx 4.1 ► CD1	•		✓ 4 Search CD1		_	
le Edit View Too	ols Help							
rganize 🔻 🛛 Burn	New folder					8== •		(
Favorites	Name	Date modified	Туре	Size				
🧮 Desktop	ISSetupPrerequisites	02/01/2018 13:44	File folder					
🐌 Downloads	🐌 Redist	02/01/2018 13:44	File folder					
🔠 Recent Places	🗿 0x0409.ini	27/07/2004 18:36	Configuration sett	6 KB				
	Autorun.inf	21/11/2005 23:29	Setup Information	1 KB				
Libraries	Data1.cab	22/11/2005 00:07	Cabinet File	483,196 KB				
Documents	instmsi30.exe	16/11/2004 06:50	Application	1,957 KB				
🌢 Music	🛃 ISScript1050.Msi	22/10/2004 03:19	Windows Installer	971 KB				
Pictures	🛱 Micromass MassLynx V4.1.msi	22/11/2005 08:58	Windows Installer	19,968 KB				
🛃 Videos	setup.exe	21/11/2005 23:29	Application	302 KB				
	Setup.ini	22/11/2005 00:07	Configuration sett	2 KB				
Computer			2					
GS (C:)								
TOSHIBA EXT (E:)								
TOSHIBA (F:)								
ms (\\128.40.177.1	.9							
Network								

There will be some questions to answer please answer as below:-

MassLynx Security Options Please select a security mode	Water
Options No Security MassLynx Basic Security MassLynx Full Security	Waters Laboratory Informatics
Install to: C:WassLynx	

😸 MassLynx V4.1 - Wizard	
MassLynx Options MassLynx Installation directory	<b>W</b> aters
Information Install MassLynx V4.1 to: C:\MassLynx	Change
InstallShield	< Back Next > Cancel
MassLynx Options Do you have a Waters 996 PDA or W	Vaters CapLC PDA system Installed?
Options Yes No	Waters Laboratory Informatics MaSSLynt
Install to: C:\MassLynx	
TestellChield	

Just click Next

Application Mana Please select the r	ager Options equired Application Mana	agers from the list below: Wate
Options BioLynx FractionLynx MaxEnt MicrobeLynx OLGS ProteinLynx TargetLynx QuanOptimize	ChromaLynx MarkerLynx MetaboLynx NeoLynx ProfileLynx PSD MX Transform	Waters Laboratory Informatics MassLyn
Install to: C:WassLynx stallShield		

## **Click Next**

😸 MassLynx V4.1 - Wizard	X
Ready to Install the Program The wizard is ready to begin installation.	Waters
Click Install to begin the installation.	
If you want to review or change any of your installation settings, click Back. C exit the wizard.	lick Cancel to
InstallShield	Cancel

### **Click Install**

HassLynx V4.1 - Wizard	
Ready to Install the Program The wizard is ready to begin installation.	<b>W</b> aters
Click Install to begin the installation.	
If you want to review or change any of your installation settings, click Back. exit the wizard.	Click Cancel to
InstallShield	
< Back Install	Cancel

If you have any questions please email to mass spec staff.