Laboratory Risk Assessment

Procedure	Removal of Carbonates Using 10% Hydrochloric Acid
-----------	---

Level of Risk	Medium
	1110 01101111

Hazard	Risk	k Recommended Control		
Preparation of 10% Hydrochloric Acid from Concentrated Acid	Personal injury - chemical burns, eye injury	?? Lab coat, safety glasses and gloves must be worn at all times?? Conc. HCl must be used in the fume cupboard with the window pulled down to the safe working height as indicated		
Addition of Acid to Samples in Test Tubes	High reactivity causing sample to "bubble over"	 ?? Always add acid to water. ?? Procedure must be carried out in a fume cupboard. ?? Always wear PPE ?? Always start by adding the acid dropwise. ?? If sample reacts violently, add water from a squeazy bottle to dilute the acid and slow the reaction down. ?? If sample spills from tube, ensure the area is washed down and all residues cleaned from the outside of the tube. 		

Assessment for C.O.S.H.H.

Procedure	Preparation of samples for diatom analysis
11000000	Treparation of Samples for diatom analysis

Substance/Procedure	Risk of exposure L/M/H*	HSE Exposure Limits (mg/m³)	Local controls used	Disposal	Emergency procedures
Preparation of 10% Hydrochloric Acid from Concentrated Acid	L	7	F/C, PPE, DG,	B, G	1, 5, 6

^{*} Risk of exposure (Low, Medium or High) providing local controls are used For key to symbols see separate table